



Using the Master Thermometer with an Auto Dialer

After attaching the power supply to the Auto Dialer:

1. Program the Master Thermometer (see Care & Use guide), to collect and display temperatures.
2. Program the Master Thermometer to trigger the auto dialer using the Master Thermometer menus
 - a. Set the alarm sound time (at least 1 minute), alarm limits and delay times on each sensor setup menu..
3. Program the auto dialer (see AVD-45 Installation & Instruction Manual) with phone numbers to call. **Set auto dialer for N/O, No exit delay, Entry delay and Mom.**
4. Attach the RJ11 phone plug from the dialer to the green relay of the Master Thermometer (white and yellow wires go into the green relay. It does not matter which wire goes where.
5. Plug an analog phone line RH11 plug into the auto dialer

Make sure that the dialer is programmed for a N/O (normally open) relay or the dialer will start calling numbers as soon as you plug it into a phone line.